

anti-ssDNA/APOSTAIN

Monoclonal Antibody. For research use only. Not for in vitro diagnostic or therapeutic procedures.

Cat.No.:BMS156

PRODUCT INFORMATION AND MANUAL



*Bender MedSystems GmbH
Campus Vienna Biocenter 2
A-1030 Vienna, Austria *Europe*

TABLE OF CONTENTS

1.	PRINCIPLES OF THE PROCEDURE	3
2.	MONOCLONAL ANTIBODY F7-26	5
3.	APPLICATIONS	6
4.	PROTOCOLS	11
5.	REFERENCES	16
6.	ORDERING INFORMATION	19

DETECTION OF APOPTOTIC CELLS WITH MONOCLONAL ANTIBODY TO SINGLE-STRANDED DNA

1. PRINCIPLES OF THE PROCEDURE

Apoptosis assay described herein is based on the increased sensitivity of DNA in apoptotic cells to thermal denaturation. In this method DNA is denatured by heating in the presence of formamide and stained with monoclonal antibody (MAb) F7-26 specific to single-stranded DNA (ssDNA).

The protocols presented below are based on the modification of the previously described procedure for the detection of apoptotic cells with MAb specific to ssDNA (1-3). The main part of this modification is the heating cells and tissue sections in the presence of formamide, instead of PBS-MgCl₂ solution (4,5).

The formamide-MAb procedure has the following advantages: 1. Selective denaturation of DNA in apoptotic cells after the heating at a relatively low temperature (56-75° C instead of 100° C).

2. Specific staining of apoptotic cells in thin sections of formalin-fixed tissues prepared with routine histological techniques (previously described procedure produced best results after methanol fixation). Indeed, this new protocol produced best and stable results after formalin fixation.

This procedure specifically stains condensed chromatin of apoptotic cells. Since chromatin condensation is a definite marker of apoptosis (6), the histochemical identification of condensed apoptotic chromatin by anti-ssDNA MAb makes possible specific detection of apoptotic cells. The specificity of the assay is based on the high sensitivity of DNA in condensed chromatin of the apoptotic cells to thermal denaturation (1,7). In the presence of formamide, only in apoptotic nuclei DNA denatures and binds the anti-ssDNA MAb F7-26 after heating at 56-75° C (4,5).

Importantly, formamide induces DNA denaturation in the apoptotic cells, and did not affect the stability of DNA with single- or double-stranded DNA breaks in the absence of apoptosis (4,5). As a result, formamide-MAb procedure specifically stains apoptotic cells and clearly distinguishes apoptosis from necrosis. Independence of DNA breaks is critical, because techniques that detect double-stranded DNA breaks, such as TUNEL, are not specific for

apoptosis (8-16). Indeed, necrotic cells in three models of necrosis were brightly stained by TUNEL, but were completely negative by formamide-MAb staining (4,5).

The following facts demonstrate that selective denaturation of DNA in apoptotic nuclei in the presence of formamide is not related to DNA breaks:

1. Necrotic cells with a high level of double-stranded DNA breaks, demonstrated by TUNEL, do not react with the anti-ssDNA MAb.

2. Apoptotic cells with very low and high levels of DNA fragmentation are stained by formamide-MAb technique with a similar intensity. 3. Nuclei with single-stranded DNA breaks induced by oxygen radicals do not bind the MAb after the heating in formamide, although alkaline denaturation detects DNA breaks in these nuclei (4,5).

These data indicate that DNA denaturation in apoptotic cells in the presence of formamide is induced by the changes in chromatin associated with the condensation per se independently of DNA breaks. As histones are known to stabilize DNA against thermal denaturation (17), it is possible that changes in DNA-histone interactions in condensed chromatin of the apoptotic nuclei may be responsible for the increased DNA sensitivity to thermal denaturation. Indeed, digestion of histones occurs during apoptosis (18,19) and increased immunoreactivity and extractability of histones precedes appearance of DNA breaks (20).

Monoclonal antibody F7-26 used in this procedure was generated against calf thymus ssDNA and selected on the basis of reactivity with mouse apoptotic cells. This antibody specifically reacts with deoxycytidine and requires for its binding ssDNA of at least 25-30 bases in length (2,21). MAb F7-26 is **not species-specific** and could be used **to detect apoptotic cells in various species (mouse, rat and human cells were tested)**.

2. MONOCLONAL ANTIBODY F7-26, BMS156

A. Catalog number

Mab F7-26 holds the catalog number BMS156.

B. Specificity

Mab F7-26 was generated by the immunisation of mice with calf thymus ssDNA and was initially used as a tool for the detection of drug-induced DNA breaks (24,25). The antibody was applied for the staining of apoptotic cells when the conditions for the selective denaturation of apoptotic DNA were developed (22,26). The most important characteristic of F7-26 is the binding to ssDNA and the absence of any reactivity with DNA in double-stranded conformation (24,25). The specific binding to ssDNA plays a critical role in the application of Mab F7-26. The anti-ssDNA Mabs F7-26 and AP-13 produce identical staining of apoptotic cells, although the antigenic determinants of these antibodies are different (deoxycytidine and deoxythymidine, respectively) (23,25).

The assay is highly specific for apoptotic cells as demonstrated by the absence of MAb reactivity with necrotic cells or cells with DNA breaks in the absence of apoptosis. Specificity of the assay was demonstrated using necrosis induced by hyperthermia or detergents and DNA breaks induced by hydrogen peroxide. These cells were completely negative by formamide-MAb assay, but positive by TUNEL (1,2). Specificity is a major advantage of anti-ssDNA MAb assay in comparison with widely used TUNEL procedure.

C. Dilution and storage


The Mab F7-26, BMS156, is supplied as 1 ml of frozen solution containing approximately 100 µg/ml of the antibody (IgM isotype) in PBS supplemented with 5% fetal bovine serum (FBS). To prepare working concentration of the Mab, thaw the ampoule in room temperature water and add 9 ml of PBS supplemented with 5% FBS. This will provide 10 ml of Mab solution sufficient for 100 tests (100 µl/test). The diluted antibody should be aliquoted and stored frozen. The antibody is stable at -20°C for 2 months and at -70°C/-80°C for 3 years. The working solution could be frozen/thawed 3 - 4 times without the loss of activity.

3. APPLICATIONS

A. FLUORESCENCE MICROSCOPY.

In cultures of breast cancer MB-MDA-468 cells treated with protein kinase inhibitor staurosporine, only cells with chromatin condensation typical of apoptosis were labeled with antibody to ssDNA. In contrast cells with dispersed chromatin in staurosporine-treated cultures or interphase nuclei of untreated cells did not react with the MAb (Fig 1A,B).

A. DNA fluorochrome DAPI

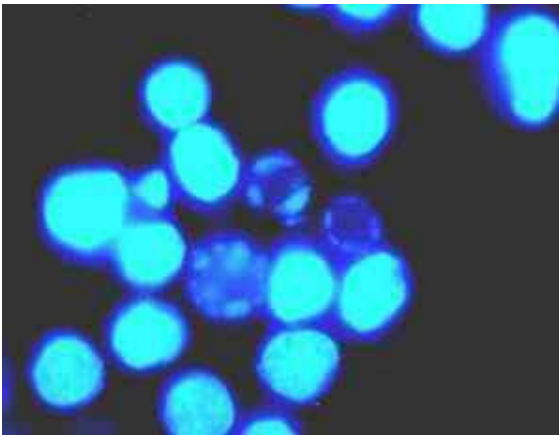

B. MAb to ssDNA

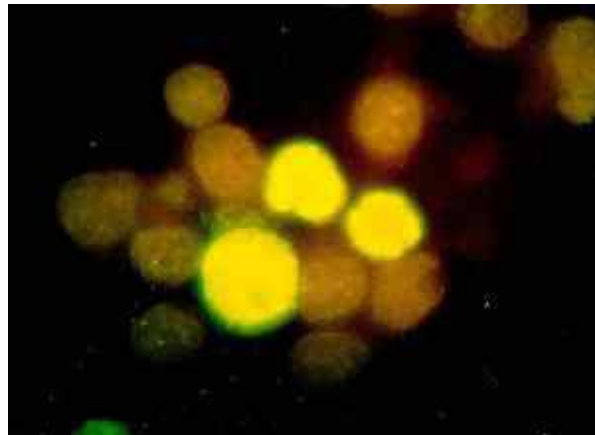


FIGURE 1. Fluorescence microphotographs of MB-MDA-468 cells treated with staurosporine, heated in formamide, stained with MAb F7-26 and counterstained with DNA fluorochrome DAPI. The same field is shown after UV excitation for DAPI (left panel) and visible light excitation for fluorescein-labeled antibody (right panels). Note that only **3 apoptotic cells** with chromatin condensed at the nuclear periphery are stained with the MAb. Magnification, x1000

B. FLOW CYTOMETRY.

The ability of formamide-MAb technique to distinguish between apoptosis and necrosis is demonstrated by flow cytometric analysis of cells treated with apoptosis and necrosis inducing agents. Fluorescence profiles of cell heated in the presence of formamide and stained with MAb to ssDNA are demonstrated in Fig. 2A. Two subsets of cells were observed in cultures treated with staurosporine: cells with low fluorescence similar to that of the untreated cells and brightly fluorescent apoptotic cells. In contrast to the high fluorescence intensity of apoptotic cells, the necrotic cells in cultures

treated with the inhibitor of respiratory chain sodium azide or with hyperthermia had fluorescence profile similar to that of untreated control cells. In contrast, TUNEL stained not only apoptotic, but also necrotic cells (see Fig 2B).

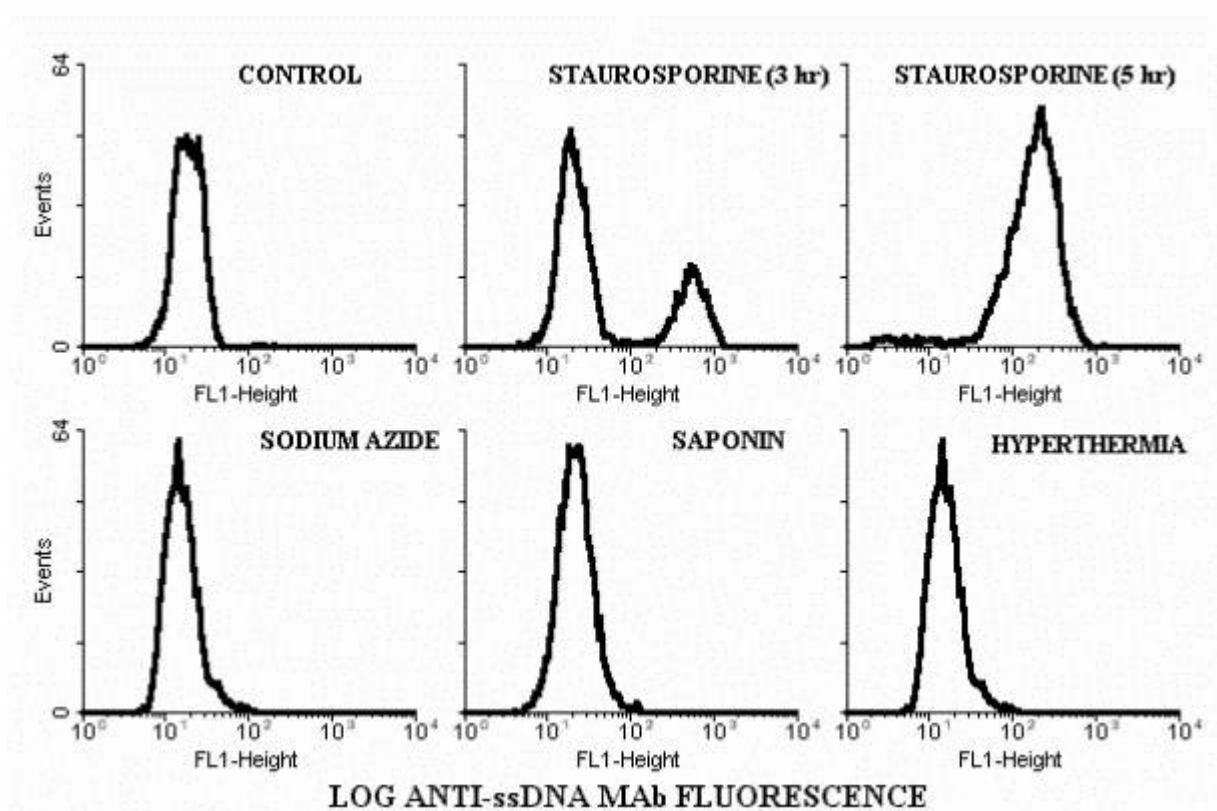


FIGURE 2A. Fluorescence distributions of MB-MDA-468 cells heated in formamide and stained with MAb F7-26 generated on a flow cytometer. Apoptosis was induced by staurosporine and necrosis was induced by sodium azide or hyperthermia. Note that apoptotic cells are intensely stained with the MAb, while fluorescence profiles of necrotic and control cells are similar.


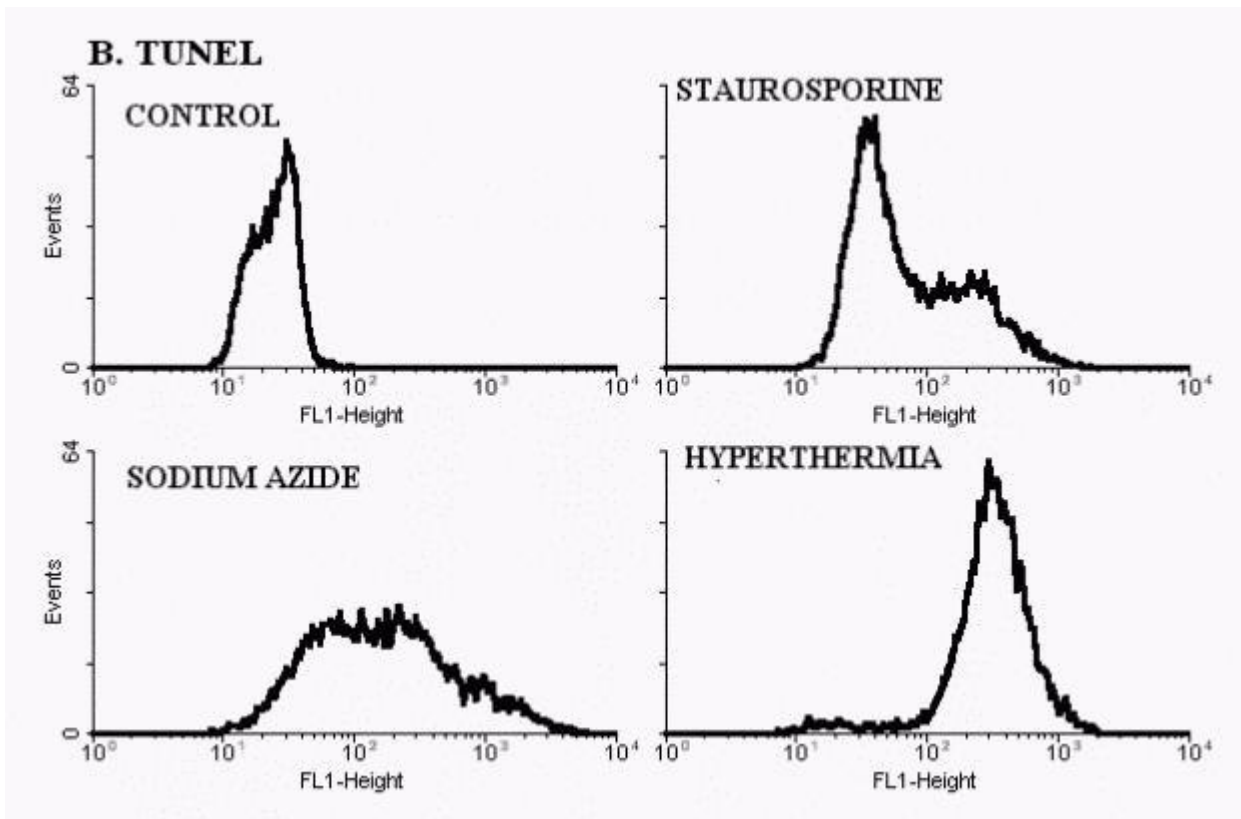


FIGURE 2B. Fluorescence distributions of MB-MDA-468 cells stained with TUNEL generated on a flow cytometer. Apoptosis was induced by staurosporine and necrosis was induced by sodium azide or hyperthermia. Note that apoptotic and necrotic cells are stained.


C. DETECTION OF APOPTOSIS IN TISSUE SECTIONS.

Small intestine of mice treated with DNA synthesis inhibitor, hydroxyurea and thymus of mice treated with hydrocortisone were fixed in formalin. Thin sections of tissues prepared with standard histological techniques were heated in formamide and stained with MAb F7-26.

Apoptotic cells with condensed chromatin and fragmented nuclei in the crypts of the small intestine from hydroxyurea-treated mice were stained with the MAb (Fig. 3A). There was no staining of cells in the crypts of untreated animals (Fig.3B). Since crypts contain large number of S-phase and mitotic cells these data indicate that this MAb does not stain normal proliferating cells.

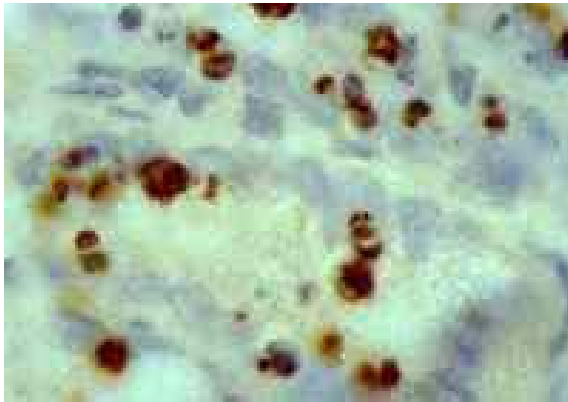

A. Hydroxyurea

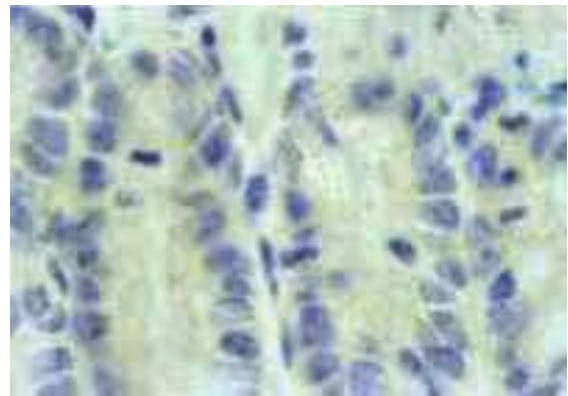


B. Control

FIGURE 3. Photomicrographs of small intestine crypts of a mouse treated with hydroxyurea (**A**) and small intestine crypts of a control animal (**B**), treated. Sections of formalin-fixed tissues were heated in formamide, stained with MAb F7-26 and counterstained with hematoxylin. Brown nuclei - antibody stained apoptotic cells with condensed chromatin and fragmented nuclei.

Multiple foci of MAb positive apoptotic cells were detected in the thymus of mice treated with the glucocorticoid, while only rare stained cells were seen in the thymus of untreated animals (Fig.4A,B).

A. Hydrocortisone

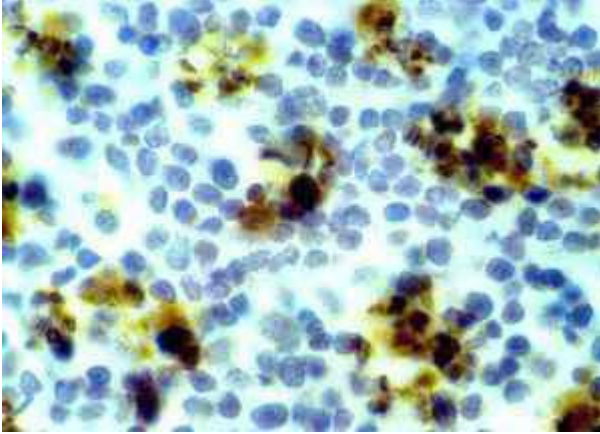

B. Control

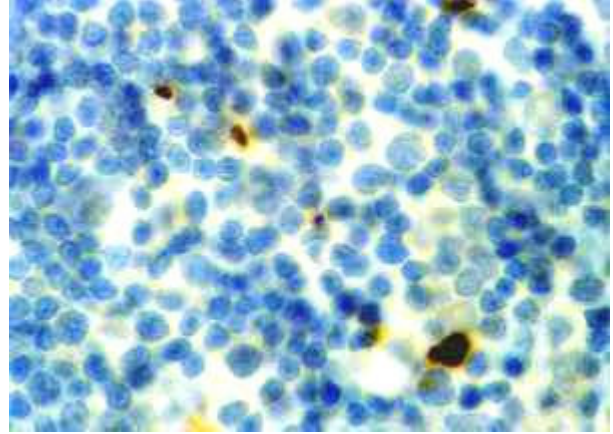


FIGURE 4. Photomicrographs of thymus of a mouse treated with hydrocortisone (**A**) and thymus of a control animal (**B**). Sections of formalin-fixed tissues were heated in formamide, stained with MAb F7-26 and counterstained with hematoxylin. Brown nuclei - antibody stained apoptotic cells with condensed chromatin and fragmented nuclei.

4. PROTOCOLS

A. Staining of tissue sections

I. Paraffin sections

Procedure includes the following steps:

1) Preparation of thin sections from **formalin** –fixed tissues using standard histological techniques. 2) Permeabilization in **saponin** and treatment with **Proteinase K**. 3) Heating in **formamide**. 4) Staining with anti-ssDNA monoclonal **antibody F7-26** and peroxidase conjugated **anti-mouse IgM**.

1. Fix tissues in 4% or 10% neutral buffered formalin (or 4% paraformaldehyde in PBS) at +4° C for 24 hours, dehydrate and embed in paraffin. **Important: fix in cold formalin.**
2. Cut 3-4 micron sections from paraffin blocks and attach sections to superfrost/plus slides. Stain freshly cut sections.
3. Deparaffinize tissue sections in xylene, wash sequentially in 100%, 95%, 70% ethanols and PBS.
4. Incubate slides in PBS containing 0.2 mg/ml saponin and 20 µg/ml Proteinase K at room temperature for 20 min.
5. Wash slides in distilled water
6. Transfer slides into coplin jar containing 40 ml of 50% formamide (v/v distilled H₂O) preheated in water bath to 56-60° C. Incubate slides in coplin jar with formamide for 20 min in water bath. **Important: temperature of formamide solution inside the jar should be 56° C.** *Formamide (Sigma Chemicals, cat.# F-7503).*
7. After the heating, transfer slides into ice-cold PBS for 5 min.
8. Quench endogenous peroxidase in 3% hydrogen peroxide for 5 min. Rinse in PBS.

9. Treat sections with 3% non-fat dry milk (Carnation) for 20 min at 37 C to block non-specific antibody binding. Rinse in PBS. *To prepare blocking solution dissolve non-fat dry milk powder in distilled water and preheat to 37 C.*
10. Apply 100 µl of monoclonal antibody F7-26 to the slide. Incubate at room temperature for 30 min and rinse in PBS. *To prepare working concentration of the antibody, add 9 ml of 1% non-fat dry milk in PBS to the vial containing 100 µg of F7-26. The diluted antibody should be aliquoted and stored frozen at -20 C.*
11. Apply 100 µl of peroxidase-conjugated anti-mouse IgM, incubate 30 min and rinse with PBS. *To prepare working concentration of the antibody, dilute peroxidase-conjugated rat monoclonal anti-mouse IgM (Zymed, cat # 04-6820) 1:100 in PBS. Use only **freshly prepared** solution. We recommend peroxidase-conjugated anti-IgM because biotin-avidin systems may produce non-specific cytoplasmic staining in some tissues. Signal intensity was strong without biotin-avidin amplification. Sigma peroxidase-conjugated goat anti-mouse IgM (cat.#A8786) also could be used as a secondary antibody.*
12. Apply chromagen solution (DAB), counterstain with hematoxylin and mount.

II. Frozen Sections

Two protocols are recommended for frozen sections

- I.
 - 1) Fix tissue in **cold** 4% paraformaldehyde in PBS for 18-24 hours.
 - 2) Cut frozen sections.
 - 3) Immerse slides with sections into 80% methanol in PBS for 30 min
 - 4) Air-dry fixed sections and store dried at room temperature.

- 5) Immerse dried sections in saponin /Proteinase K solution as in step 4 for paraffin sections and proceed as in protocol described above.
- II.**
- 1) Cut frozen sections from fresh tissue
 - 2) Fix sections in 4% paraformaldehyde in PBS for 10 min
 - 3) Rinse in PBS
 - 4) Immerse slides with sections into 80% methanol in PBS for 30 min
 - 5) Air-dry fixed sections and store dried at room temperature.
 - 6) Incubate slides with dried sections in PBS containing 0.2 mg/ml saponin for 20 min (Proteinase K is not used in this protocol)
 - 7) Proceed to step 5 for paraffin sections.

B. Staining of archival formalin-fixed tissues

The standard protocol described above provides optimal results after fixation in cold formalin or paraformaldehyde. Fixation at room temperature or prolong fixation may inhibit staining, because excessive DNA-protein cross-links prevent formamide-induced DNA denaturation in apoptotic cells.

Additional protease (Pronase E) treatment is recommended for archival material if weak or negative staining is obtained with the standard protocol.

1. After Saponin/Proteinase K treatment rinse slides in PBS and Tris buffer (10 mM, pH 7.6)
2. Incubate slides in Pronase E (20 microg/ml Tris buffer) at room temperature for 20 min in Coplin jar. *Pronase E (Sigma Chemicals, cat.#P6911) is prepared as a stock solution 1mg/ml Tris buffer and stored frozen.*

3. Rinse slides in distilled water and proceed to the heating in formamide as in standard protocol.

C. Staining of cell suspensions for flow cytometry and fluorescence microscopy

Procedure includes the following steps

- 1) Fixation in **methanol**.
- 2) Heating in **formamide**.
- 3) Staining with anti-ssDNA monoclonal **antibody F7-26** and fluorescein-conjugated **anti-mouse IgM**.
 1. Resuspend cell pellets in 1 ml of cold PBS and slowly add 6 ml of methanol precooled to -20°C while vortexing.
 2. Store fixed cell at -15 - 20°C for 1-3 days before staining.
 3. Transfer 5×10^5 of fixed cells into 15 ml plastic tubes (Falcon 2096), centrifuge, remove supernatant.
 4. Resuspend pellet in 0.25 ml of formamide. Keep 5 min at room temperature.
 5. Immerse rack with tubes into water bath preheated to 75 degrees C for 10 min. *Circulating water baths are recommended for optimal heating.*
 6. After the heating, immediately transfer rack with tubes into room temperature water
 7. 2 ml of 1% non-fat dry milk in PBS to the tubes containing 100% formamide, vortex and keep tubes at room temperature for 15 min.
 8. Centrifuge, remove supernatant, resuspend pellet in 100 microL of monoclonal antibody F7-26 and incubate at room temperature for 15 min. *To prepare working concentration of the antibody, add 9 ml of 5% fetal*

bovine serum in PBS to the vial containing 100 micro g of F7-26. The diluted antibody could be aliquoted and stored frozen at -20 °C.

9. Add 1 ml of PBS, centrifuge and resuspend pellet in 100 microL of fluorescein-conjugated anti-mouse IgM. Incubate 15 min at room temperature. *To prepare working concentration of the antibody, dilute fluorescein-cojugated goat anti-mouse IgM (Sigma, Cat. # F-9259) in 1% non-fat dry milk in PBS. The diluted antibody could be aliquoted and stored frozen at -20 °C.*
10. For flow cytometry, add 1 ml of PBS, centrifuge and resuspend pellet in 0.5 ml of propidium iodide solution (1 microg/ml in PBS).
11. For fluorescence microscopy, rinse cells with PBS, prepare cytopspins and stain slides with DNA fluorochrome DAPI (0.1 microg/ml in PBS for 10 min).

8. REFERENCES

1. Frankfurt, O.S. (1994) Detection of apoptotic leukemic and breast cancer cells with monoclonal antibody to single-stranded DNA. *Anticancer Res.* 14:1861-1870.
2. Frankfurt, O.S., Robb J. A., Sugarbaker E. V. and Villa L. (1996) Monoclonal antibody to single-stranded DNA is a specific and sensitive cellular marker of apoptosis *Exp. Cell Res.* 226:387-397.
3. Frankfurt, O. S., Robb, J.A., Sugarbaker E.V., and Villa, L. (1997) Apoptosis in breast carcinomas detected with monoclonal antibody to single-stranded DNA: Relation to bcl-2 expression, hormone receptors and lymph node metastases. *Clin. Cancer Res.* 3:465-471.
4. Frankfurt, O.S. and Krishan, A. (2001) Identification of apoptotic cells by formamide induced DNA denaturation in condensed chromatin. *J. Histochem. Cytochem.* 49:369-378.
5. Frankfurt, O.S. and Krishan, A. (2001) Enzyme-linked immunosorbent assay (ELISA) for the specific detection of apoptotic cells and its application for rapid drug screening. *J. Immunol. Methods.* 253: 133-144.
6. Zamzani, N., and Kroemer, J. (1999) Condensed matter in cell death. *Nature* 401:127-128.
7. Allera, C., Lazzarini, G., Patrone, E., Alberti, I., Barboro, P., Sanna P., Melchiori, A., Parodi, S., and Balbi, C. (1997) The condensation of chromatin in apoptotic thymocytes shows a specific structural change. *J. Biol. Chem.* 272:10817-10822
8. Grasl-Kraipp, B., Ruttkau-Nedecky, B., Koudelka, H., Bukowska, K., Bursch, W., and Schulte-Hermann, R. (1995) *In Situ* detection of fragmented DNA (TUNEL assay) fails to discriminate among apoptosis, necrosis and autolytic cell death: A cautionary note. *Hepatology* 21:1465-1468.
9. Charriaut-Marlangue, C., and Ben-Ari, J. (1995) A cautionary note on the use of the TUNEL stain to determine apoptosis. *Neuroreport* 7:61-64
10. Stadelmann C., Bruck, W., Bancher, C., Jellinger, K., and Lassmann, H., (1998) Alzheimer disease: DNA fragmentation indicates increased

neuronal vulnerability, but not apoptosis. *J. Neurophatol. Exp Neurol.* 57:456-464.

11. Didenko, V. V., and Hornsby P. J. (1996) Presence of double-stranded breaks with single-base 3' overhang sin cells undergoing apoptosis but not necrosis. *J. Cell Biol.* 135:1369-1376.
12. Ohno, M., Takemura, G., Ohno, A., Misao, J., Hayakawa, Y., Minatoguchi, S Fujiwara, T., and Fujiwara, H. (1998). *Circulation* 98: 1442-1430.
13. Collins, R.J., Harmon B. V., Gobe G. C., and Kerr J. F. R. (1992) Internucleosomal DNA cleavage should not be the sole criterion for identifying apoptosis. *Int. J. Rad. Biol.* 61:451-453.
14. Fukuda, K., KojiroM., and Chiu, J.F. (1993) Demonstration of extensive chromatin cleavage in transplanted Morris hepatoma 7777 tissue: apoptosis or necrosis? *Amer. J. Pathol.* 142: 935-946.
15. Dong, Z., Saikumar, P., Wenberg, J. M., and Venkatachalam, M. A. (1997) Internucleosomal DNA cleavage triggered by plasma membrane damage during necrotic cell death. *Amer. J. Pathol.* 151: 1205-1213.
16. Kanoh, M., Takemura, G., Misao, J., Aoyama, T., Nishigaki, K., and Noda, T. (1999) Significance of myocytes with positive in situ end labeling (TUNEL) in hearts with dilated cardiomyopathy. *Circulation* 99:2757-2764.
17. Tsai, Y. h., Ansevin, A. T., and Hnilica, L. S. (1975) Association of tissue specific histones with deoxyribonucleic acid *Biochemistry* 14: 1257-1264.
18. Kaufmann, S. H. (1989) Induction of endonucleolytic DNA cleavage in human leukemia cells by etoposide, camptotecin and other cytotoxic anticancer agents: a cautionary note. *Cancer Res.* 49:5870-5878.
19. Kuribayashi, N., Sakagami, H., and Takeda, M. (1996) Chromatin structure and endonuclease sensitivity in human leukemic cell lines. *Anticancer Res.* 16:1225-1230.
20. Zunino, S. J., Singh, M. K., Bass, J., and Picker, L. J. (1996) Immunodetection of histone epitopes correlates with early stages of apoptosis in activated human peripheral T lymphocytes. *Amer. J. Pathol.* 149:653-663.

21. Frankfurt, O.S. (1990) Decreased stability of DNA in cells treated with alkylating agents. *Exp. Cell Res.* 191:181-185.
22. Desjardins, L.M. and J.P. MacManus. (1995) An adherent cell model to study different stages of apoptosis. *Exp. Cell Res.* 216:380-387.
23. Eichhorn, G.L. (1962) Metal ions as stabilizers and destabilizers of deoxyribonucleic acid structure. *Nature* 194:474-475.
24. Frankfurt, O.S. (1987) Detection of DNA damage in individual cells by flow cytometric analysis using anti-ssDNA monoclonal antibody. *Exp. Cell Res.* 170:369-380.
25. Frankfurt, O.S. (1990) Decreased DNA stability in the cells treated with alkylating agents. *Exp. Cell Res.* 191:181-185.
26. Frankfurt, O.S., J.J. Byrnes, D. Seckinger, and E.V. Sugarbaker. (1993) Apoptosis (programmed cell death) and the evaluation of chemosensitivity in chronic lymphocytic leukemia and lymphoma. *Oncol. Res.* 5:37-42

9. ORDERING INFORMATION

For technical information and orders please contact:

Bender MedSystems GmbH
Campus Vienna Biocenter 2
A-1030 Vienna, Austria*Europe
phone: +43 1 796 40 40 ext. 100
fax: +43 1 796 40 40 ext. 400

Bender MedSystems, Inc.
849 Hinckley Road
Burlingame, CA 94010, USA
toll-free phone: +1 (866) 952 2112
toll-free fax: +1 (877) 952 2112
phone: +1 (650) 952 2112
fax: +1 (650) 952 2252

e-mail: techserv@bendermedsystems.com
www.bendermedsystems.com

Cat. No. BMS156 anti-ssDNA/APOSTAIN